



Tobii Pro Glasses 2

Quick Start Guide

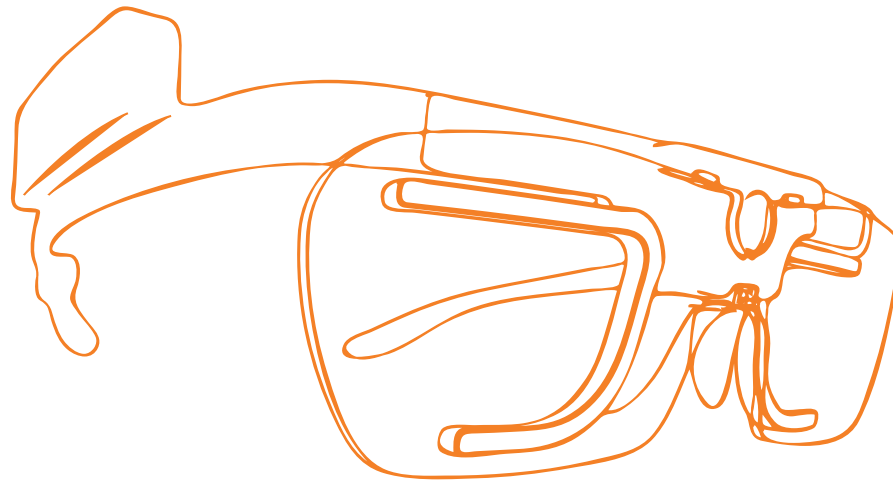

What's in the Carry Case?

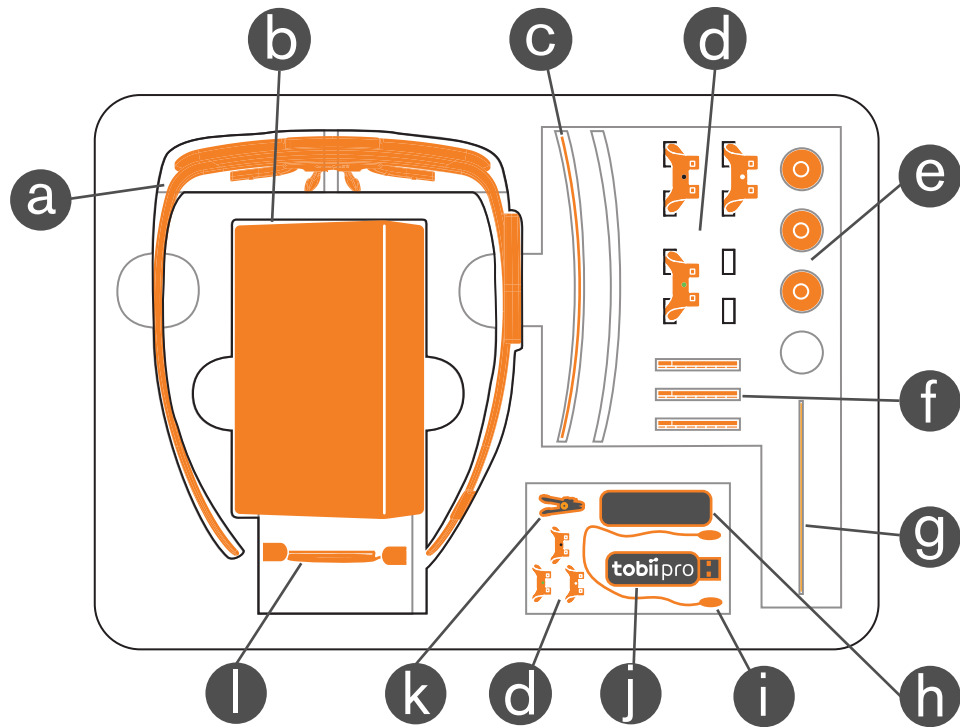

- a** 1 x Head Unit
- b** 1 x Recording Unit
- c** 2 x Protective lenses (1 clear + 1 tinted; 1 attached to the Head Unit)
- d** 6 x Nose Pads in 3 different sizes
- e** 3 x Rechargeable Li-Ion batteries type 18650
- f** 3 x SD memory cards + sleeves
- g** 1 x Calibration card
- h** 1 x Microfiber cleaning cloth
- i** 1 x Head Strap
- j** 1 x USB memory stick with documentation and software
- k** 1 x Cable cord clip
- l** 1 x HDMI A to HDMI D cable

What's in the Accessories Box?




- 1** 1 x Battery charger for Li-Ion batteries
- 2** 1 x USB Ethernet Adapter
- 3** 5 x Power cables for battery charger (EU, US, UK + AUS)
- 4** 1 x USB charger for Recording Unit
- 5** 4 x Power plugs for USB charger (EU, US, UK + AUS)
- 6** 1 x Ethernet cable (3 Meters)
- 7** 2 x Calibration cards
- 8** 10 x Calibration stickers

Preparing and Connecting Tobii Pro Glasses 2


1 - Charge the batteries

Insert the batteries into the charger. The battery is fully charged when the icon below is displayed on the screen:

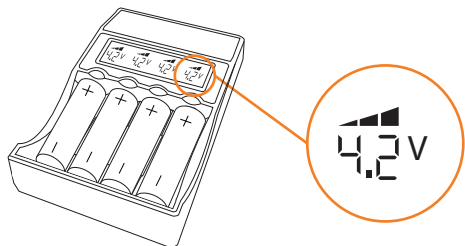

2 - Recording Unit preparation


Insert SD memory card.

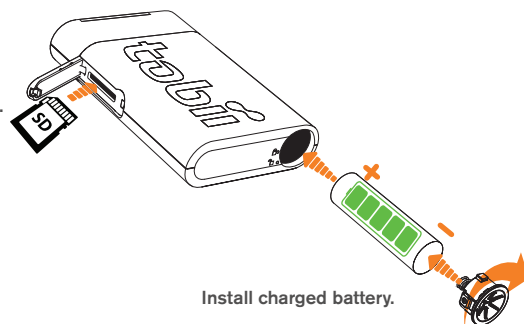

Install charged battery.

3 - Attach the Nose Pad


Gently squeeze the nose pads together and insert the Nose Pad into the two holes on the head unit.

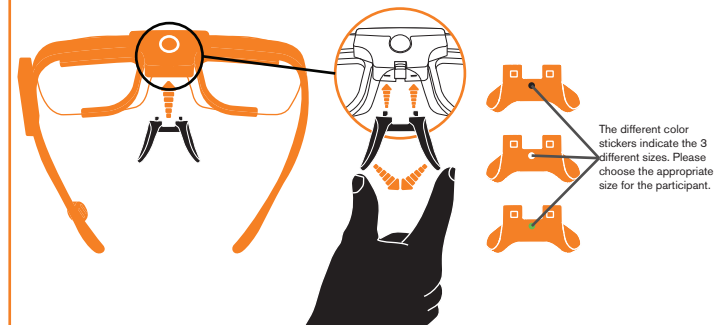


4 - Head Unit preparation

Step 1

Insert HDMI cable.

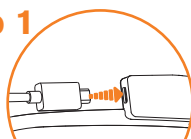


Step 2

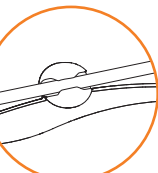

Secure cable in clip.

Optional Head Strap.

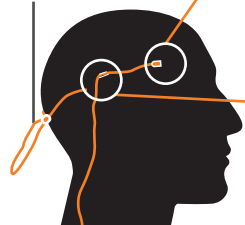


5 - Powering the Recording Unit



Insert the HDMI cable, then press and hold the power button for a few seconds or until the LEDs stop flashing to turn on your Recording Unit.

6 - Software installation

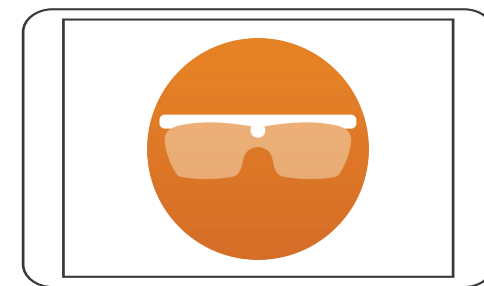

Download, install and open "Tobii Pro Glasses Controller" on your Windows tablet/computer.

www.tobii.com/downloads

Wireless

Find and connect to your Pro Glasses 2 via WLAN. (Look on the back of the Recording Unit for the serial number, this will be the name of the WLAN network.)

Step 1



The WLAN password for your Tobii Pro Glasses 2 is "TobiiGlasses" with a capital T and G.



Step 2



7 - Connecting to your device

Wired

If using a wired connection, securely connect the Ethernet cable to the Recording Unit and to your computer.

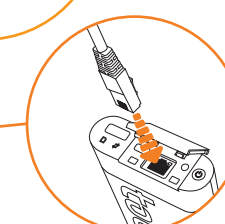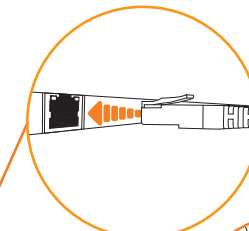

Recording Unit LEDs

	(SD LED off) The SD Memory Card ready or the Recording Unit is powered off.		(Connection LED solid green) The Recording Unit is connected to Tobii Pro Glasses Controller.
	(SD LED blinking green) The Recording Unit is busy writing to the SD Memory Card. Do NOT remove the SD Memory Card.		(Connection LED off) The Recording Unit is disconnected from Tobii Pro Glasses Controller.
	(SD LED blinking red) Missing SD Memory Card. Insert an SD Memory Card.		(Power LED off) The Recording Unit is turned off. All other LED indicators are off.
	(SD LED solid yellow) The SD Memory Card has critically low storage. SD Memory Card change recommended.		(Power LED solid green) The battery is good or charging.
	(SD LED blinking yellow) The Recording Unit is busy writing to an SD Memory Card which has critically low storage. Stop recording and change the SD Memory Card.		(Power LED solid yellow) The battery running out of power. Battery change recommended
	(SD LED solid red) The SD Memory Card is corrupted. Change the SD memory card.		(LED blinking yellow) The battery is critical. Change the battery.
	(SD LED blinking red) The SD Memory Card is Read-Only. Remove lock or replace the SD Memory Card.		(LED Slow blinking) No power to the Head Unit. Ensure the head unit is connected correctly.

1 - Creating a new recording

Start up the "Tobii Pro Glasses Controller". If this is your first ever recording, create a new Project before proceeding. Then press "Start New Recording" on the relevant pair of Tobii Pro Glasses 2.

Type the participant's name and relevant info, then press "Create".

1 - Software installation

Step 1
Download, install and open "Tobii Pro Lab Analyzer Edition" on your Windows tablet/computer.

Step 2
Remove the SD memory card from your Recording Unit and insert it into your computer.

www.tobiipro.com/downloads

2 - Calibrating a participant

Step 1
Gently fit the Head Unit onto the participant as shown below. Exchange the Nose Pad if necessary.

0.50m - 1.50m (1.65ft - 4.9ft)

Step 2
Gently attach the Recording Unit to the participant's belt or clothing.

Step 3
Hold the calibration card and have the participant focus on the center point..

Step 4
...then press the calibrate button and allow the calibration process to complete. You are now ready to start recording.

3 - Start a recording

If the calibration is successful, press the record button.


(If not, ask the participant to look at the calibration target again then press the recalibrate button.)

4 - Stop a recording

To stop a recording, press the stop button.

You have now completed a recording with the Tobii Pro Glasses 2.

2 - Create a project

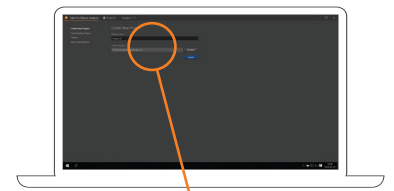

Analyze | ▾


Create New Project

Project name
Project X

Folder location
C:\Users\sse\Desktop\Project X

Using the "Tobii Pro Lab Analyzer Edition" choose a name for your new project and where to save it, then press create.

3 - Import your data

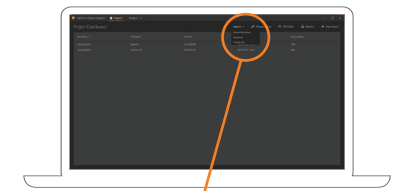

Import ▾

- Glasses Recording
- Snapshots
- Coding data

At the Project Dashboard press "Import > Glasses Recording". Locate and select the data file (*.ttgp) on the SD memory card to import it.

You are now ready to begin analyzing your data.

Please see the USB memory stick included in your package for software, a full User Manual, and an instructional video. These are also available online at:

www.tobii.com/support

Tobii_QuickStart_TobiiProGlasses2QuickStartGuide_24042017_ENGus_WEB



©Tobii®. Illustrations and specifications do not necessarily apply to products and services offered in each local market. Technical specifications are subject to change without prior notice. All other trademarks are the property of their respective owners.

Tobii Pro Support

EUROPE / GLOBAL
Phone (SWE): +46 8 522 950 10
Phone (GER): +49 69 24 75 03 4-27
support@tobii.com
Support hours: 8 am - 6 pm
Between July-August: 9am - 5pm
(Central European Time, GMT +1)

NORTH AMERICA
Phone: +1 703 738 1320
support.us@tobii.com
Support hours: 8 am - 8 pm
(US Eastern Standard Time, GMT -6)

JAPAN
Phone: +81 3 5793 3316
support.jp@tobii.com
Support hours: 9 am - 5.30 pm
(Japan Standard Time, GMT +9)

CHINA
Phone: +86 180 1558 5168
support.cn@tobii.com